
Sanitary Sensor 4113s

Pharmaceutical sanitary open type for ultrapure injection water

4-electrode conductivity sensor with external electrodes. Used with TCU series 3100 and 3300.

MaterialsPTFE, silicone rubber and steel 1.4404 (AISI 316L)Wet SS parts of sensor are pol- ished according to the pharma- ceutical standard, Ra < 0.80 μmPressure12 bar at 150 °C1.0 ± 0.5 % (electrically adjuste Calibrated in a Ø51 mm SS pipe	ı. d)
ished according to the pharma- ceutical standard, Ra < 0.80 μπ Pressure 12 bar at 150 °C 1.0 ± 0.5 % (electrically adjuste	d)
1.0 ± 0.5 % (electrically adjuste	1
	1
Calibrated in a (751 mm 55 pipe	
Cell constant Ø63 mm SS pipe gives an error of ~ -6 %.	
Ø63 mm plastic pipe gives an error of ~ 20–30 %.	
Reproduceability ±2 %	
Linearity From 20000 to 0.2 µS/cm ± 2 % per decade.	Ď
Below 5 μS/cm + 0.01 μS/cm.	
Temp. sensor Pt 1000 (IEC 751 class A)	
Time delay T ₉₀ , 9 sec.	
Protection IP65 splash-proof (DIN 40 050)	
Sensor cable PVC 12 x 0.25 mm ²	
standard length 3 m	
max. temp. 70 °C	
Sensor can optionally be mount ed with 5 m of cable; measuring below 0.5 μS/cm with this cable length is not recommended.	1
Connection 11-pole strip connector with MF20 cable adapter	
Weight 0.6 kg	

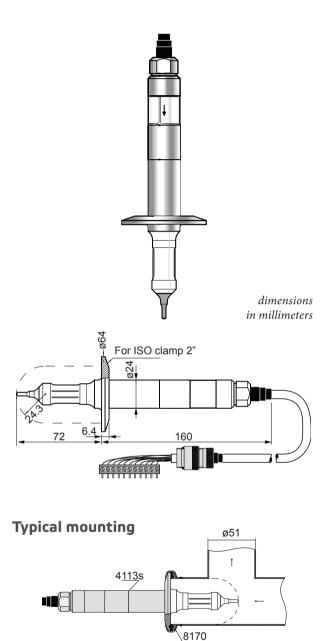

8951

4<u>113</u>s


Mounting


Sensor mounting with a 2" ISO 2852 clamp (not included). Keep a minimum distance of 24.3 mm to the surroundings.

On a horizontal pipe the sensor should be mounted to the side of the pipe, in a place where no air bubbles occur. Construct the piping in a self-draining manner.




Mounting with type 8151 stainless steel nipple in a 2" plastic Tee

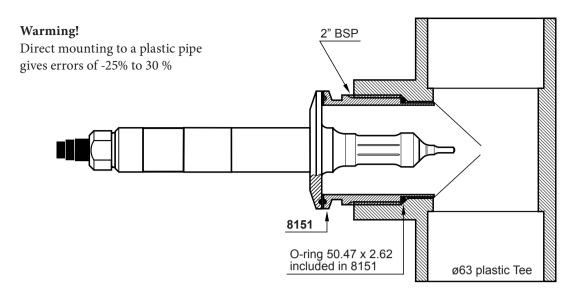

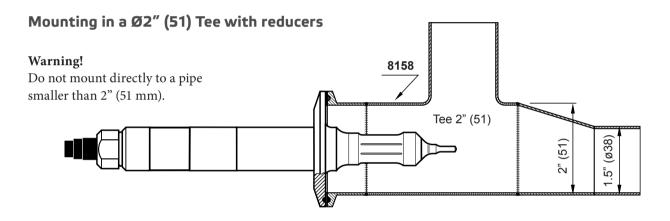

Mounting with 8155 stainless nipple in a 63.5mm stainless Tee

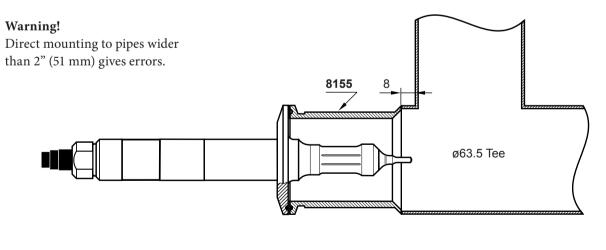

For more information, contact your local Valmet office. www.valmet.com



Specifications in this document are subject to change without notice. Product names in this publication are all trademarks of Valmet Corporation.